TIARIS™ Tissue Protect Buffer

(Molecular Biology Grade)

Ordering Info

TBB0430. TIARIS™ Tissue Protect Buffer, 50 x 1.5 mL

TBB0432. TIARIS™ Tissue Protect Buffer, 500 mL

TBB0431. TIARIS™ Tissue Protect Buffer, 100 mL

Description

TIARIS™ Tissue Protect Buffer is an effective nucleic acids stabilization and storage solution. It preserves high-quantity and good-quality RNA and DNA. Once the sample is introduced into TIARIS™ Tissue Protect Buffer, RNases are immediately inactivated so nucleic acids are preserved for long periods until the samples can be processed.

Storage

Store at 25°C.

☞ If any precipitate is observed, warm the buffer at 37°C.
Agitate to redissolve it.

Quality Control

Buffer composition and pH are verified from lot to lot.

Features

- Non-toxic solution, and easy to transport to collect field samples.
- Tissues can be stored indefinitely at -20°C or -80°C.
- Suitable for most tissues (ear, mouse tail tip, organs, etc) cultured cells, bacteria, yeast, plants.
- Compatible with most RNA isolation systems.

Applications

- RNA/ DNA Purification.
- PCR/ RT-PCR.
- Library preparation.
- Northern blot.
- Nuclease protection assay.

PROTOCOL

; Wear gloves during sample preparation !

1. Prepare the samples:

Tissue: Cut a small piece of fresh tissue (~ 0.5 cm) and cut in smaller pieces.

Cell Cultured: Pellet cells (~ 5 x 10⁶ cells) at 200 g for 5 minutes at room temperature. Wash the cells in PBS 1x pH=7.4.

Unicellular microorganisms: Pellet cells (~ 3 x 10⁸ cells) at 12,000 g for 2 minutes.

2. Add 5-10 volumes for tissue / 1 mL for cell pellets of TIARIS™ Tissue Protect Buffer.
3. Incubate the samples overnight at 4°C and store the samples attending at:

Temperature	Storage	Recommendations
-80 °C	Indefinitely	Centrifugated at 12,000 g, 5 min and discard the supernatant before store at -80°C.
-20 °C	Indefinitely	
4 °C	1 month	
25 °C	1 week	
37 °C	24 hours	

TIARIS™ Tissue Protect Buffer

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Referencias

TBB0430. TIARIS™ Tissue Protect Buffer, 50 x 1,5 mL

TBB0432. TIARIS™ Tissue Protect Buffer, 500 mL

TBB0431. TIARIS™ Tissue Protect Buffer, 100 mL

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PROTOCOLO

¡ Wear gloves during sample preparation !

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